




# DAM Series Rackmount

Model 3484

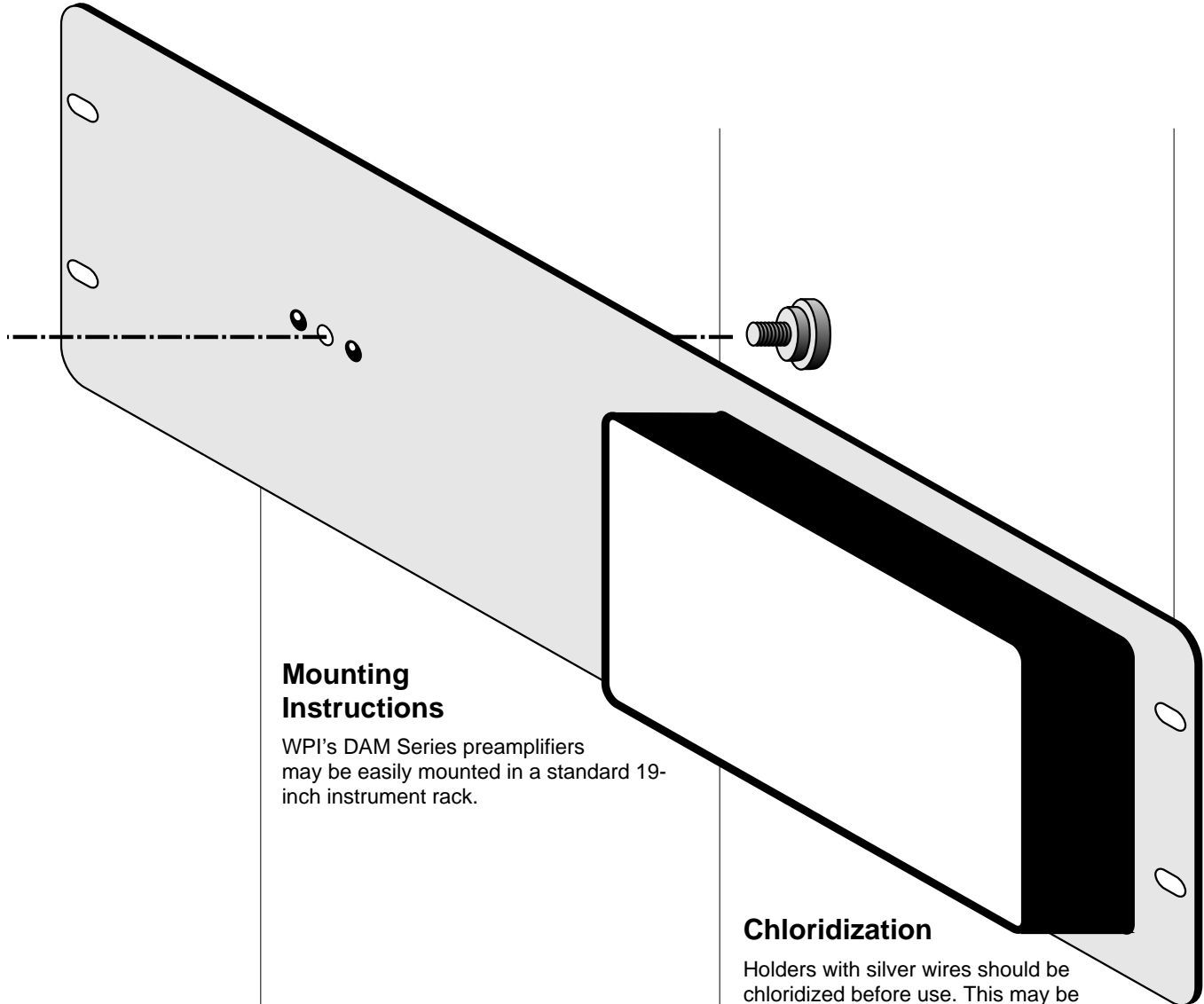

## Mounting Instructions

WPI's DAM Series preamplifiers may be easily mounted in a standard 19-inch instrument rack.

## Chloridization

Holders with silver wires should be chloridized before use. This may be done easily by immersing the wire in 1 M hydrochloric acid and applying +1.5 V to the holder vs. a return electrode, such as a large platinum wire. If, before chloridizing, the wire is tarnished (yellow to dark red) it may be cleaned by passing it quickly over a flame, being careful to avoid melting the acrylic body, or washing in urea solution.

10/91

## ***World Precision Instruments, Inc.***

International Trade Center, 175 Sarasota Center Boulevard, Sarasota FL 34240-9258 USA

Telephone: 941-371-1003 • Fax: 941-377-5428 • E-mail: [wpi@wpiinc.com](mailto:wpi@wpiinc.com)

**UK:** Astonbury Farm Business Centre, Astonbury, Stevenage, Hertfordshire SG2 7EG England • Tel: 0438-880025 • Fax: 0438-880026

**Germany:** Liegnitzer Str. 15, D-10999 Berlin, Germany • Tel: 030-6188845 • Fax: 030-6188670

**Japan:** 1-4-2 Naka-Meguro, Meguro, Tokyo 153, Japan • Tel: 03-3220-5200 • Fax: 03-3220-5201