Boom Stand

Instruction Manual



Please read the Instruction Manual carefully before installation and keep it for future use.

Open the package and check the parts

- 1. Base
- 2. Vertical post
- 3. Base locking screw
- 4. Guard ring
- 5. Cross block
- 6. Horizontal bar
- 7. Groove locking screw
- 8. Top cap
- 9. Locking screw

Please contact the local dealer for any missing part or problem.



Installation

Step 1		Step 2	Step 3
Use the Screw 3, to attach the Base 1 and the Vertical		Put the Guard Ring 4 on	Put the Cross Block 5 on
Post 2 together, and put it on the table.		the vertical post.	the Vertical post, and
			tighten the right side
			locking screw.
3-2		4	5

Step 4	Step 5
Put the Horizontal Bar 6 into the Cross Block, and tighten the left side of locking	Put the locking screw 9 on
screw. Confirm the slot of the Horizontal Bar should be on the top side, and use the	the side of Horizontal Bar.
screw 7 to fix the Horizontal Bar. When the horizontal bar is on the end of cross block,	
it can't be pulled out.	
the grown is the top	

Step 6	The End		
Put the Cap 8 on the top of the	Your installation work is done		
vertical post.	successfully!		
8			

Attention

Please check your ordered stand in the following models based on your purchase order.



Boom Stand Accessories

The following parts are not included in the boom stand package. Please contact the local dealer for the parts.

76mm/83mm / 85mm	39mm	Fine focus E-Arm (for any	B&L
E-Arm	E-Arm	Dia. of the scope holder)	E-Arm
-	-		

CCD Mounting Plate	CCD Mounting Plate with	76mm Through Hole	32mm / 25mm Arbor
	Focus	Focus Rack	
-			Ť

How to attach the Standard 76mm E-Arm			
On the top	On the front of	On the bottom. Warning: please use the safety screw	
	Horizontal Bar	when you put the E-Arm on the bottom of the bar.	
			-