

# 1

## UNBOX

The scope head and stand are shipped separately. Carefully unbox the microscope head.

**TIP:** Firmly grasp the poly bag to pull the scope head out of the box. DO NOT LIFT the head by the visor.



# 2

## STAND

Assemble the stand according to the manufacturer's directions.

**NOTE:** The stand and scope are shipped in separate boxes.



**ESTIMATED  
SETUP TIME:  
2 min.**

# 3

## MOUNT SCOPE

Line up the focus mount bracket on the back of the head with the bracket on the stand.

Slide the head into place.


The Universal Quick Connect comes with your stand.



# 4

## POWER

Plug the power cord into the wall outlet, and plug the other end of the cable into the power port on the back of the head.

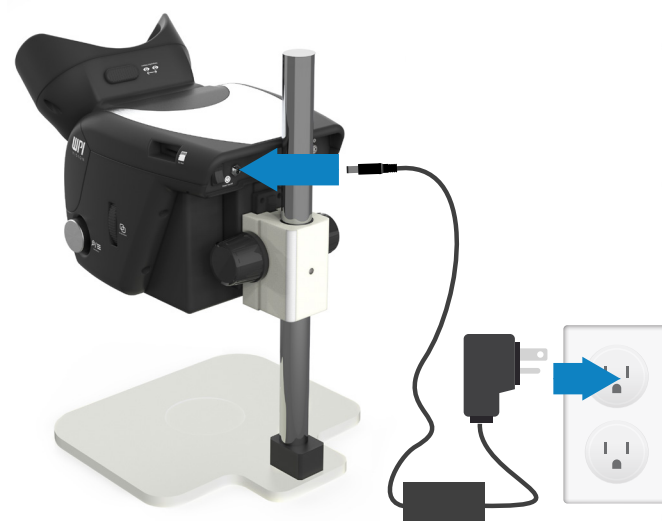

# 5

## FUNCTION KNOBS & BUTTONS

- Press and release either one of the Image buttons to capture a still image. Press and hold an Image button to take video. Release the button to stop the video capture.
- Press either Menu/Zoom button to access the on-screen menu. Rotate the knob to zoom.

- Rotate the Light Intensity knob to adjust the lighting.

The Menu/Zoom and Image buttons are redundant for ambidextrous use.

Light Intensity  
Menu/Zoom  
Picture  
Menu/Zoom  
Video



# 6

## ALL SET

You are ready to make your work life easier, faster, more comfortable and accurate. Enjoy!

**TIP:** To view your saved pictures and videos on a personal device we recommend using VLC software.



# SAVE THE BOX.

Please save the packaging, in case you need to return the unit.

A Division of

**WORLD PRECISION INSTRUMENTS**

USA: 175 Sarasota Center Boulevard, Sarasota FL 34240-9258 USA • Tel: (941) 371-1003 • Fax: (941) 377-5428 • wpi@wpiinc.com • www.wpiinc.com

UK: 1 Hunting Gate, Hitchin, Hertfordshire SG4 0TJ England • 44 (0)1462 424700 • wpiuk@wpi-europe.com

Germany: Pfingstweide 16, D-61169 Friedberg (Hessen) • Tel: +49 (0)6031 1602171 • Fax: +49 (0)6031 1602180 • wpide@wpi-europe.com

China and Hong Kong: Rm 27A, No8 Donfang Rd., Pudong District, Shanghai 200120 PRC • +86 688 85517 • ChinaSales@china.wpiinc.com