



HS6

Micromanipulator

CARE & MAINTENANCE

This micromanipulator is a precision built instrument. It was calibrated at the factory to standard parameters and is ready for use. As with any delicate mechanical device, it needs your best care and attention to ensure many years of accurate and trouble-free performance.

Please handle this instrument with care. Dropping it even a short distance can be damaging to the general alignment and adjustment. Dust is a major source of problems for this instrument. When not in use, adjust the three (X-Y-Z) guide surfaces in such a way that they are not exposed and cover the manipulator with a plastic bag. If not in use for a prolonged time, occasionally move the three (X-Y-Z) surfaces in a series of forward and reverse motions to keep the grease pliable. Never oil the guide surfaces. Under normal use, you should not need to apply grease for many years.



World Precision Instruments, Inc.



WARRANTY

WPI (World Precision Instruments, Inc.) warrants to the original purchaser that this equipment, including its components and parts, shall be free from defects in material and workmanship for a period of one year* from the date of receipt. WPI's obligation under this warranty shall be limited to repair or replacement, at WPI's option, of the equipment or defective components or parts upon receipt thereof f.o.b. WPI, Sarasota, Florida U.S.A. Return of a repaired instrument shall be f.o.b. Sarasota.

The above warranty is contingent upon normal usage and does not cover products which have been modified without WPI's approval or which have been subjected to unusual physical or electrical stress or on which the original identification marks have been removed or altered. The above warranty will not apply if adjustment, repair or parts replacement is required because of accident, neglect, misuse, failure of electric power, air conditioning, humidity control, or causes other than normal and ordinary usage.

To the extent that any of its equipment is furnished by a manufacturer other than WPI, the foregoing warranty shall be applicable only to the extent of the warranty furnished by such other manufacturer. This warranty will not apply to appearance terms, such as knobs, handles, dials or the like.

The foregoing obligations set forth in this paragraph are in lieu of all obligations and liabilities, including all warranties of merchantability or otherwise, expressed or implied or statutory, and state WPI's entire and exclusive liability and purchaser's exclusive remedy for any claim of damages in connection with the sale or furnishing of all equipment, including design, suitability for use, operation, or installation. There are no warranties which extend beyond the description of the face hereof. In no event shall WPI be liable for any special or consequential damages.

Warning: This equipment is not designed or intended for use on humans.

** Electrodes, batteries and other consumable parts are warranted for 30 days only from the date on which the customer receives these items.*

World Precision Instruments, Inc.

USA: International Trade Center, 175 Sarasota Center Boulevard, Sarasota FL 34240-9258 USA
Tel: 941-371-1003 • Fax: 941-377-5428 • E-mail: sales@wpiinc.com • Internet: <http://www.wpiinc.com>

UK: Astonbury Farm Business Centre • Aston, Stevenage, Hertfordshire SG2 7EG England • Tel: 01438-880025 • Fax: 01438-880026 • E-mail: wpiuk@wpi-europe.com

Germany: Liegnitzer Str. 15, D-10999 Berlin, Germany • Tel: 030-6188845 • Fax: 030-6188670 • E-mail: wpide@wpi-europe.com

Japan: 3-6-14-202 Kasumigaseki, Chiyoda, Tokyo 100-0013, Japan • Tel: 81-3-3592-4500 • Fax: 81-3-3592-4540 • E-mail: sales@wpi-j.com

Australia: P.O. Box 1191, Glen Waverley, Victoria 3150, Australia • Tel: (03) 9887-6262 • Fax: (03) 9887-9585 • E-mail: wpi.au@ozemail.com.au