

24-Hour Emergency Number (ChemTel) 888.255.3924 in U.S.

cleaning@alconox.com

alconox.com

- distributors
- technical information
- free samples
- new developments
- white papers
- prices



ALCONOX® Powdered Precision Cleaner

- Concentrated to save you money
- Biodegradable and readily disposable
- Penetrating wetting power to save you time
- Replaces corrosive acids and hazardous solvents
- Free rinsing to provide reliable results and no interfering residues
- Use to pass cleaning validation tests for lab accreditation and plant inspection approval

Used to clean: Healthcare instruments, laboratory ware, vacuum equipment, tissue culture ware, personal protective equipment, sampling apparatus, catheters, tubing, pipes, radioactive contaminated articles, optical parts, electronic components, pharmaceutical apparatus, cosmetics manufacturing equipment, metal castings, forgings and stampings, industrial parts, tanks and reactors. Authorized by USDA for use in federally inspected meat and poultry plants. Passes inhibitory residue test for water analysis. FDA certified.

Used to remove: Soil, grit, grime, buffing compound, slime, grease, oils, blood, tissue, salts, deposits, particulates, solvents, chemicals, radioisotopes, radioactive contaminations, silicon oils, mold release agents.

Surfaces cleaned: Corrosion inhibited formulation recommended for glass, metal, stainless steel, porcelain, ceramic, plastic, rubber and fiberglass. Can be used on soft metals such as copper, aluminum, zinc and magnesium if rinsed promptly. Corrosion testing may be advisable.

Cleaning method: Soak, brush, sponge, cloth, ultrasonic, flow through clean-in-place. Will foam—not for spray or machine use.

Directions: Make a fresh 1% solution (2 1/2 tbsp per gal, 1 1/4 oz per gal, or 10 g per L) in cold, warm, or hot water. If available, use warm water. Use cold water for blood stains. For difficult soils, raise water temperature and use more detergent. Clean by soak, circulate, wipe, or ultrasonic method. Not for spray machines: will foam.

For nonabrasive scouring, make paste. Use 2% solution to soak frozen stopcocks. To remove silver tarnish, soak in 1% solution in aluminum container. RINSE THOROUGHLY—preferably with running water. For critical cleaning, do final or all rinsing in distilled, deionized, or purified water. For food contact surfaces, rinse with potable water. Used on a wide range of glass, ceramic, plastic, and metal surfaces. Corrosion testing may be advisable. Material compatibility guidelines available on request.

Convenient Sizes	Alconox Cat. #	
Case 9 x 4 lb (9 x 1.8 kg) boxes	1104	A
25 lb (11 kg) carton	1125	
50 lb (22 kg) carton	1150	
100 lb (45 kg) drum	1101	
300 lb (136 kg) drum	1103	
Case 12 box 50 x 1/2 oz (15 g)	1112	No. 1
1 lb makes 13 gal (1 kg makes 1 cleaning solution	00L)	

ALCONOX is available from leading laboratory, hospital, clinical and industrial suppliers. To find a distributor for Alconox, Inc. detergents, use the Locate Dealers search tool in the Support section of alconox. com. To request FREE samples, use the Get Samples form at alconox.com. Or call Alconox, Inc. at +914.948.4040, today.

AXTB.1





24-Hour Emergency Number (ChemTel) 888.255.3924 in U.S.

cleaning@alconox.com

alconox.com

- distributors
- technical information
- free samples
- new developments
- white papers
- prices

Physical Data		Typical Value
pH of 1% Solution		9.5
Surface Tension 1% Sol'n (dyne/cm)		32
Conductivity 1% Sol'n	(dyne/cm at 21 degrees C) 7.0
Form		Powder
Color	White and cream colored flakes	
Solubility in Water	To 10% (w/w) at ambient temperature	
Hard Water Effectiven	ess	Highly effective
Biodegradability	Yes, mee	ets EC648/2004
Foam Tendency		High foaming
Shelf Life 24	months from end of month	of manufacture
Fragrance Content		0%
Phosphate Content (as	phosphorus)	7.3%
Flash Point (degrees F)		None
Organic Carbon (% calc	ulated w/w)	11%

Chemical Description

ALCONOX consists primarily of a homogeneous blend of sodium linear alkylaryl sulfonate, alcohol sulfate, phosphates, carbonates. ALCONOX is anionic in nature.

Cleaning Validation Methods

To quantify trace levels of Alconox for cleaning validation purposes, use conductivity, UV, TOC, HPLC, or phosphorus detection. For details see the Cleaning Validation page in Support at alconox.com. For lab accreditation tests to show the detergent is rinsed away, see the Accreditation Residue Detection in the Support section of alconox.com.

Health Safety Information

OSHA Hazardous Ingredie	ents None	
RCRA Hazard Class	Nonhazardous	
EPA Priority Pollutants	None	
DOT Hazard Class	Nonhazardous	
Flammability	Nonflammable	
Latex Content	None in product or packaging.	
Irritation	Cat. 2 skin, Cat. 2A eye irritant	
Oral Toxicity	$LD_{50} > 5000 \text{ mg/kg oral rat.}$	
Inhalation Toxicity	Solution: No, Powder dust: irritant	
VOC Content	0%	
GMO Content	None	
BSE/TSE Source	None	
2007/68/EC Annex IIIa Allergens		
REACH Art. 57 (SVHC)	None	
RoHS	No RoHS hazardous ingredients.	
Carcinogenicity	NTP = No IARC = No OSHA = No	
Indgredients on TSCA Inv	rentory Yes	
EC 907/2006 Co	ontains no enzymes, optical brighteners,	
perfumes, allerg	jenic fragrances, or preservative agents.	

Precautions

No special precautions other than good industrial hygiene and safety practices employed with any industrial chemical (see Directions). A Safety Data Sheet is available at alconox.com.

Contact Alconox, Inc. for purchase specifications. Information presented is typical. Not to be taken as specifications. Typical data is not a specification.

While the information in this report should not be considered to be a product warranty, we urge you to investigate, test and verify the suitability of Alconox, Inc. detergents for your specific application. We, of course, can not give permission to use, or recommend the use of, our detergents where they infringe patents. No representation or warranty is made as to the safety of products or materials mentioned under the Federal Food Additives Amendment of 1958.