

### Warranty

WPI (World Precision Instruments, Inc.) warrants to the original purchaser that this equipment, including its components and parts, shall be free from defects in material and workmanship for a period of 6 months from the date of receipt. WPI's obligation under this warranty shall be limited to repair or replacement, at WPI's option, of the equipment or defective components or parts upon receipt thereof f.o.b. WPI, Sarasota, Florida U.S.A. Return of a repaired instrument shall be f.o.b. Sarasota.

The above warranty is contingent upon normal usage and does not cover products which have been modified without WPI's approval or which have been subjected to unusual physical or electrical stress or on which the original identification marks have been removed or altered. The above warranty will not apply if adjustment, repair or parts replacement is required because of accident, neglect, misuse, failure of electric power, air conditioning, humidity control, or causes other than normal and ordinary usage.

To the extent that any of its equipment is furnished by a manufacturer other than WPI, the foregoing warranty shall be applicable only to the extent of the warranty furnished by such other manufacturer. This warranty will not apply to appearance terms, such as knobs, handles, dials or the like.

WPI makes no warranty of any kind, express or implied or statutory, including without limitation any warranties of merchantability and/or fitness for a particular purpose. WPI shall not be liable for any damages, whether direct, indirect, special or consequential arising from a failure of this product to operate in the manner desired by the user. WPI shall not be liable for any damage to data or property that may be caused directly or indirectly by use of this product.

### Claims and Returns

- Inspect all shipments upon receipt. Missing cartons or obvious damage to cartons should be noted on the delivery receipt before signing. Concealed loss or damage should be reported at once to the carrier and an inspection requested. All claims for shortage or damage must be made within 10 days after receipt of shipment. Claims for lost shipments must be made within 30 days of invoice or other notification of shipment.

Please save damaged or pilfered cartons until claim settles. In some instances, photographic documentation may be required. Some items are time sensitive; WPI assumes no extended warranty or any liability for use beyond the date specified on the container.

- WPI cannot be held responsible for items damaged in shipment en route to us. Please enclose merchandise in its original shipping container to avoid damage from handling. We recommend that you insure merchandise when shipping. The customer is responsible for paying shipping expenses including adequate insurance on all items returned.

- Do not return any goods to WPI without obtaining prior approval and instructions (RMA#) from our returns department. Goods returned unauthorized or by collect freight may be refused. The RMA# must be clearly displayed on the outside of the box, or the package will not be accepted. Please contact the RMA department for a request form.

- Goods returned for repair must be reasonably clean and free of hazardous materials.

- A handling fee is charged for goods returned for exchange or credit. This fee may add up to 25% of the sale price depending on the condition of the item. Goods ordered in error are also subject to the handling fee.

- Equipment which was built as a special order cannot be returned.

- Always refer to the RMA# when contacting WPI to obtain a status of your returned item.

- For any other issues regarding a claim or return, please contact the RMA department

**Warning: This equipment is not designed or intended for use on humans.**

## Kwik-Cast™ & Kwik-Sil™

*Silicone elastomers cure in five minutes*

Serial No. \_\_\_\_\_



**WARNING: MAY CAUSE AN ALLERGIC SKIN REACTION.**

If the silicone elastomers make skin contact, wash with plenty of soap and water. If skin irritation or rash occurs, seek medical advice or attention. Remove contaminated clothing and wash before reuse. Wash thoroughly after handling this product. Wear protective gloves, protective clothing, eye protection and face protection.

[www.wpiinc.com](http://www.wpiinc.com)

#### USA

175 Sarasota Center Blvd., Sarasota FL 34240-9258  
Tel: 941-371-1003 • Fax: 941-377-5428 • E-mail: sales@wpiinc.com

#### UK

1 Hunting Gate, Hitchin, Hertfordshire SG4 0TJ  
Tel: 44 (0)1462 424700 • Fax: 44 (0)1462 424701 • E-mail: wpiuk@wpi-europe.com

#### Germany

Saarstraße 23, D-61169 Friedberg (Hesson), Germany  
Tel: +49 (0)6031 1602171 • Fax: +49 (0)6031 1602180 • E-mail: wpide@wpi-europe.com

#### China & Hong Kong

WPI Shanghai Trading Co., Ltd.  
Rm 18A, No8 Dongfang Rd., Pudong District, Shanghai, 200120 PR China  
Tel: +86 21 6888 5517 • E-mail: chinasaless@china.wpiinc.com

#### Brazil

Av. Conselheiro Nébias, 756 sala 2611, Santos-CEP: 11045-002, São Paulo Brazil  
Tel: (013) 406-29703 • E-mail: info@brazil.wpiinc.com

#### Internet

[www.wpiinc.com](http://www.wpiinc.com) • [www.wpi-europe.com](http://www.wpi-europe.com) • [www.wpiinc.net](http://www.wpiinc.net) • [www.wpi brasil.com.br](http://www.wpi brasil.com.br)

**Kwik-Sil** and **Kwik-Cast** silicones are based on technology with vinyl terminated siloxane and platinum complex catalysts. In order to gain enough speed to cure at room temperature, special cross linkers and high catalyst concentration is used. Although the high concentration of catalysts makes these products more costly than traditional RTV silicones and less attractive for general usage, they provide an excellent value in applications for the biological research field.



Fig. 1–Kwik-Sil and Kwik-Cast are packaged in an easy-to-use dual-syringe applicator which eliminates the introduction of air bubbles during mixing and causes a more consistent curing time.

Both **Kwik-Sil** and **Kwik-Cast** have very low toxicity before, during and after curing. In traditional RTV silicone systems, a by-product of the condensation (curing) is either acetic acid or alcohol, which are toxic to living cells. In vinyl systems, the by-product of condensation is a small amount of hydrogen gas, which is much less toxic to the cell.

**Kwik-Sil** and **Kwik-Cast** curing speed is hundreds of times faster than traditional RTV silicones. A curing time of a few minutes at room temperature is especially useful for encapsulation of live tissue or implanting into a live animal.

Unlike many vinyl-based silicones in which the platinum complex catalysts are easily poisoned by contamination from amines and animal tissue, **Kwik-Sil** and **Kwik-Cast** are not sensitive to contamination from animal tissue.

### Kwik-Sil

**Kwik-Sil** is a translucent, medium-viscosity silicone adhesive, developed for chronic peripheral nerve studies such as anterograde tracing with fluorescent indicators or electrode recording. Good adhesion and mechanical properties (tear strength and elongation) allow days of study without breaking of the bonding. Curing speed is very reproducible.

### Kwik-Cast

**Kwik-Cast** is a very low viscosity silicone sealant developed to embed peripheral nerves with electrodes for acute multi-fiber recordings. It flows easily, filling the small spaces around the nerve and leaving no channels through which peritoneal fluid can travel and thus short the nerve/electrode contact. Equally important is the ability of the material to flow into itself and create one continuous mass from underneath the nerve all the way to the top of the nerve/electrode contact to ensure long-term recording stability.

**Kwik-Cast** is color-coded to make the mixing foolproof. The catalyst is yellow and the base is blue. When uniformly mixed, it is green. **Kwik-Cast** can be applied and cured underneath mineral oil. After recording, electrodes are easily recovered due to the low tear strength.

### Storing

Store below 27°C. Keep the applicator capped to avoid air contamination when not in use.

### Mixing and Application

**NOTE:** To remove air bubbles from tip area, place capped syringe upright for several hours or over-night before use.

The cap is keyed so that it can only be applied in one direction, to avoid cross-contamination of catalyst and base. The new dual-syringe applicator mixes the silicone as you apply it.

1. Remove the white protective cap from the syringe.
2. Slip on a mixer tip (supplied in a separate bag) and turn it 90° to lock the tip in position. As the silicone is extruded from the syringe and applied directly onto the subject, it is automatically mixed by the tip. (The disposable mixer tip can only be used once because the silicone inside becomes solid very quickly. Only a very small amount of silicone is wasted in the tip, because its dead volume is less than 0.12 mL.)
3. After dispensing the silicone, the mixer tip may be left on the syringe until next application and replaced by a new one. However, if the silicone is not going to be used for several days, we recommend covering the syringe with the protective cap. Please note that the cap is keyed so that it can only be put on in one direction.

### Curing

Silicones are formulated to cure with a number of catalysts, and each catalyst cures under its own set of conditions. The curing process may be inhibited by your material preparation, device components or even assembly processes. The “poisoning” could leave a slight surface tack or cause a complete failure to cure. In some cases, the effects can be reversed, but often the inhibitors permanently arrest the cure.

The platinum catalyst used in the **Kwik-Sil** and **Kwik-Cast** silicones is susceptible to poisoning. The following list details the most common inhibitors (but not all):

#### Sulfur containing compounds

- Latex rubber (gloves & tubing)
- Neoprene rubber
- Natural Rubber
- Polysulfides

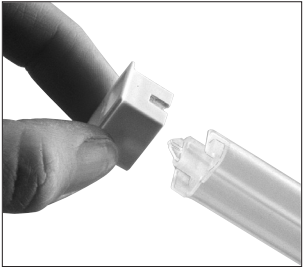


Fig. 2–Remove protective cap from syringe. After use, remove nozzle and re-cap.

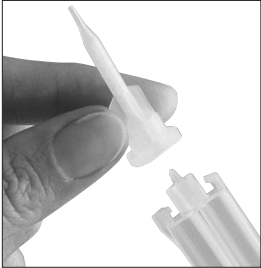


Fig. 3–Baffles in applicator nozzle ensure thorough mixing of expressed silicone.

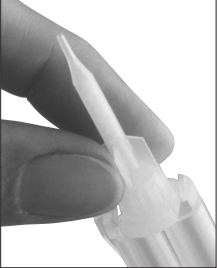


Fig. 4–Twist nozzle to lock into position.

- Wood (spatulas and tongue depressors)
- Plasticized PVC
- Organic oils

#### Epoxy & amine cured materials

- Some phenolic resins
- Epoxy resins
- Some polyurethane materials

#### Nitrogen containing compounds

- Amines
- Amides
- Nitriles
- Cyanates

#### Tin containing compounds

- Condensation cured silicones
- Acetoxy cured silicones
- Oxime cured silicones
- Tin Plasticizers (some plastics and PVC)

**Metals** - Various metals including silver, tin, lead and mercury

### Specifications

UNMIXED:	Kwik-Sil <sup>1</sup>	Kwik-Cast
Mix Ratio	1:1	1:1
Working time	<5 min. <sup>2</sup>	4 min.
Setting time (room temperature, 1:1 ratio)	5–10 min. <sup>3</sup>	<10 min.
Cure time	~15 min.	
Viscosity, cps	15,000	10,000
Shelf life at 23 °C	6 months	6 months
Volume	5 mL	5 mL
Number of mixing tip <sup>4</sup>	10	10
Dead volume of the mixing tip	<0.12 mL	<0.12 mL

#### AFTER CURING 24 HOURS:

Tear Strength, ppi	90	44
Elongation %	650	60
Durometer (shore A-2)	30	36
Color	translucent	green
Volume Resistivity	1x10 <sup>15</sup> Ω/cm <sup>2</sup>	1x10 <sup>15</sup> Ω/cm <sup>2</sup>

<sup>1</sup> 0–90 sec.	Liquid State	mixed silicone is able to flow under gravity
90–180 sec., <5 min.	Working Time	glue is firm, but still workable
5–10 min.	Firm Set	glue may be sticky
15–18 min.	Full Set	silicone is cured and is not sticky

<sup>2</sup> Three minutes is average, with about 90 seconds of liquidity.

<sup>3</sup> No longer mixable at this point.

<sup>4</sup>Ten mixer tips are included. If additional mixer tips are needed, order WPI part #600009 (package of 10).

### Sterilization

The silicone adhesive inside the syringes cannot be sterilized by any means including UV, gamma, H<sub>2</sub>O<sub>2</sub>, autoclaving or EtO. The exterior and the tips can be immersed in or wiped with 70/30 alcohol, or you can use a liquid sterilant like Cidex OPA (WPI #7364-4) or Rapicide OPA (WPI #504611).